

## Operator's Manual



## The Next Rockers

### Congratulations!

Congratulations on your purchase of a Next Rocker by Next Advance, Inc.


Please read this operator's manual which explains proper operation of the instrument. This manual can be viewed on the web at [http://www.nextadvance.com/manuals/Next\\_Rocker\\_Manual.htm](http://www.nextadvance.com/manuals/Next_Rocker_Manual.htm).

### Contents


Parts of the Next Rocker .....	2
Setup .....	2
Operation.....	2
Troubleshooting .....	2
Support.....	2
Warranty .....	2

# OPERATOR'S MANUAL FOR THE NEXT ROCKER

## PARTS OF THE NEXT ROCKER

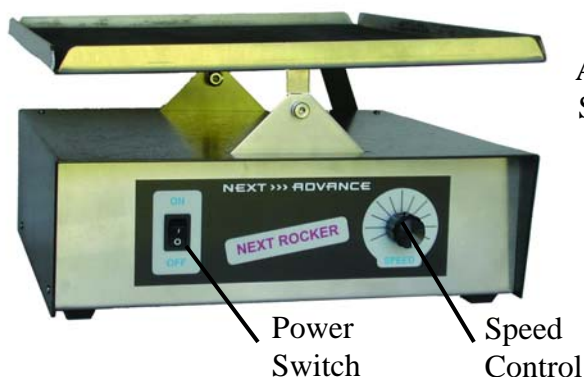

*Front of Rocker*

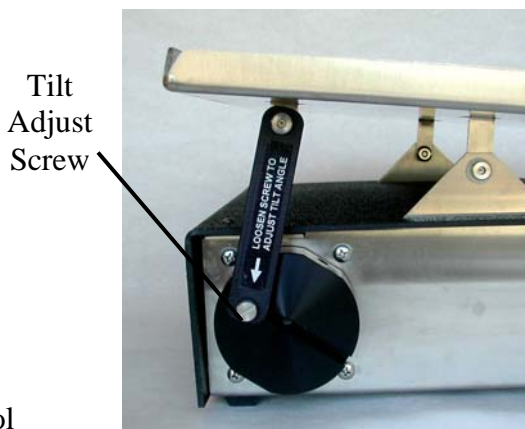

*Rear of Rocker*

## SETUP

Place the Next Rocker on a level lab bench. Plug the power supply into the back of the Next Rocker and then plug it into a wall outlet. It is now set up.

## OPERATION

Turn on the Power Switch and turn the Speed Control knob to adjust the rocking rate.

To adjust the tilt angle, use a flat blade screwdriver to loosen the labeled Tilt Adjust Screw on the rear of the instrument, slide the linkage to the desired position, then tighten the screw.

## TROUBLESHOOTING

In addition to the tips given below, a thorough list of troubleshooting tips is at [http://www.nextadvance.com/FAQs/FAQs-Next\\_Rocker.htm](http://www.nextadvance.com/FAQs/FAQs-Next_Rocker.htm).

**If nothing happens**, the power supply plug might not be in a live wall outlet or the power supply connector plug might not be fully in the jack on the back of the Next Rocker.

**If the unit stops working**, turn the system off for 15 minutes to allow the electronics to reset. If the Next Rocker does not turn on after this period, contact customer service.

## SUPPORT

A thorough list of FAQs, is at [http://www.nextadvance.com/FAQs/FAQs-Next\\_Rocker.htm](http://www.nextadvance.com/FAQs/FAQs-Next_Rocker.htm).

If you cannot find an answer there, please contact customer service by email at [support@nextadvance.com](mailto:support@nextadvance.com) or by telephone.

## WARRANTY

The Next Rocker comes with a 30 day money back guarantee (less shipping charges) and a one year warranty. Next Advance will replace, free of charge, any part which is defective due to workmanship or materials. For further information, go to [http://www.nextadvance.com/legal\\_terms.htm](http://www.nextadvance.com/legal_terms.htm). In no event shall Next Advance be liable for incidental or consequential damages.

### Warranty is Void if

Product has defect or damage due to product accident, alteration, connection to an improper electrical supply, fire, flood, lightning, or other conditions beyond the control of Next Advance.

Product is improperly installed or used.

### Operator's Responsibility

Provide proof of purchase and provide normal care and maintenance.