



QUICK INSTALLATION GUIDE FOR CRYOLYS® EVOLUTION

1. Unpacking

1. Open the box and pull out the instrument using the central handle



2. Remove the foam around the instrument and check the integrity of the unit



3. Get the unit out of the box and proceed to installation



2. Installation

Required tool: n°5 Allen key

WARNING ! Unplug the Precellys® Evolution power supply before to proceed.

1. Unscrew and remove the Precellys® Evolution lid



2. Place the new lid on the Precellys® Evolution without the cooling module



3. Insert the reinforced hinge over the Precellys® hinge and screw it to the lid



4. Plug the 2 wires at the rear of the Precellys® Evolution



5. Attach the wires to the hinge using the fast straps



6. Update the software using the USB device included



WARNING! Be careful to the number of pins of the plugs